



EQUIPMENT CALIBRATION VERIFICATION

(Ref. [SOP VAL-175](#))

Project Number	Autoclave Number
----------------	------------------

1. Procedure

Equipment calibration verification must be performed prior to the commencement of autoclave validation.

1.1 Calibration Checks

1.1.1 Verify the autoclave is within calibration (VAL-220). Print out Equipment Maintenance Status Report (Engineering Department) and attach to this document.

Sign/Date

1.1.2 Complete a list of test equipment used (Table 1). Attach a copy of calibration certificates to this document.

Sign/Date

Table 1. Test Equipment

Equipment Name	Serial Number	Calibrated (Yes/No)	Certificate Attached
Pressure Transducer			
Eurotherm Chessel 6000 Graphic Recorder			
Hart Scientific HDCR Calibrator			
Graphtec GL midi logger CL220/820			

1.1.3 Calibrate thermocouples (VAL-210). Copy Calibration Verification Record (Form 825) and attach to this document.

Sign/Date

1.1.4 Sign and attach all printouts of the thermocouple calibrations to this document.

Sign/Date

2. Completion

Upon completion of this test attach document to the Autoclave Validation Protocol (Form 780).