



AUTOCLAVE CYCLE PARAMETERS RECORDING SHEET - DRY

(Ref. [SOP VAL-175](#))

DRY - CYCLE # Parameters

Verify that all cycle parameters are set as listed in the table below.

Parameter Description		Expected Parameter Set Point	Actual Parameter Set Point	Acceptance Criteria Met (Yes/No)
Cycle Number		(Enter number)		
Cycle Name		(Enter Name)		
DRY-PreC: Evacuate 1	Value (kPa)	≤ 30		
	Delay (s)	60		
DRY-PreC: Charge 1	Value (kPa)	≥ 150		
	Delay (s)	60		
DRY-PreC: Evacuate 2	Value (kPa)	≤ 25		
	Delay (s)	60		
DRY-PreC: Charge 2	Value (kPa)	≥ 150		
	Delay (s)	30		
DRY-PreC: Evacuate 3	Value (kPa)	≤ 20		
	Delay (s)	60		
DRY-PreC: Charge 3	Value (kPa)	≥ 150		
	Delay (s)	1		
Heating Ramp-Up to S.P (pressure)	Value (kPa)	≥ 219.3kPa (equiv. 123.0°C)		
	Delay (s)	1		
DRY: Sterilisation	Value (min)	≥ 15		
	Delay (s)	1		
Chamber Exhaust	Value (kPa)	≤ 120		
	Delay (s)	1		
Vacuum Drying	Value (min)	≥ 10		
	Delay (s)	1		
Chamber Equalisation	Value (kPa)	≥ 110		
	Delay (s)	60		
Vacuum Drying	Value (min)	≥ 15		
	Delay (s)	1		
Chamber Equalisation	Value (kPa)	≥ 110		
	Delay (s)	60		
Vacuum Drying	Value (min)	≥ 10		
	Delay (s)	1		
Chamber Equalisation	Value (kPa)	≥ 105		
	Delay (s)	1		
Sterilisation Control Method	Chamber Pressure @	Chamber Pressure @ 219.3kPa		

Action	Acceptable (Yes/No)	Sign/Date	Verification Sign/Date
All cycle parameters are set correctly			